

WinCamD

The Camera we are using belongs to the Chip experiment.

Technical Data

TaperCamD20-15-UCD23:

- WinCamD with 20 x 15 mm 2.27:1 FO taper on the sensor
- 2/3" CCD sensor for CW Pulsed, 1360 x 1024 pixels, 6.45 μm

If you have a TaperCamD series camera, press **Alt S** to reach the WinCamD Capture Setup dialog, and enter the appropriate PMF value in the WinCamD capture screen.

- TaperCamD20 - 15 - UCD23 Pixel multiply factor = **2.27**

Also enter the correct wavelength.

Congratulations. You have successfully completed installation.



Shortcuts

Start measurement: Press G

Stop measurement: Press S

Power Limits

Power $> 2.5 \times (\text{Beam diam in mm}) \text{ W}$, or $> 10 \text{ W}$ total, may damage the ND filter. The limit is lower for pulsed lasers, especially ns & fs lasers.

nm	Damage Threshold	Saturation Irradiance
	mW/cm ²	uW/cm ²
355	Few	~0.3
630	>10	~0.15
800	"	~0.1
1060	"	~10

Data sheet

- WinCamD: [wincamd_man.pdf](#)

Last
update:
2015/11/30 13:51 groups:mg:experiment:cameras:wincam <https://iqwiki.iqo.uni-hannover.de/doku.php?id=groups:mg:experiment:cameras:wincam&rev=1448891460>

From:
<https://iqwiki.iqo.uni-hannover.de/> - IQwiki

Permanent link:
<https://iqwiki.iqo.uni-hannover.de/doku.php?id=groups:mg:experiment:cameras:wincam&rev=1448891460>

Last update: **2015/11/30 13:51**

