



Fiberlaser

Close the program "*Fiber Comb Control*" and lock out the CEO- & REP-PID in the rack, otherwise you can't optimize the comb!

- Push Yes

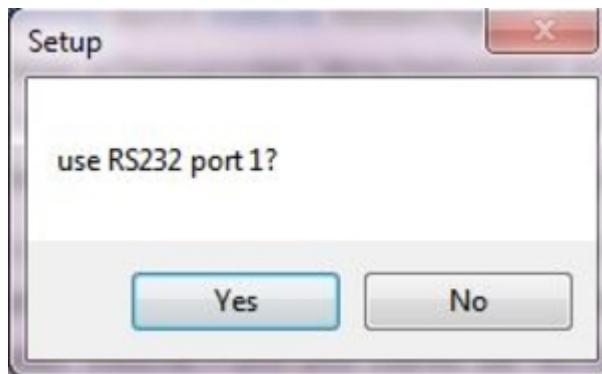

- Push OK

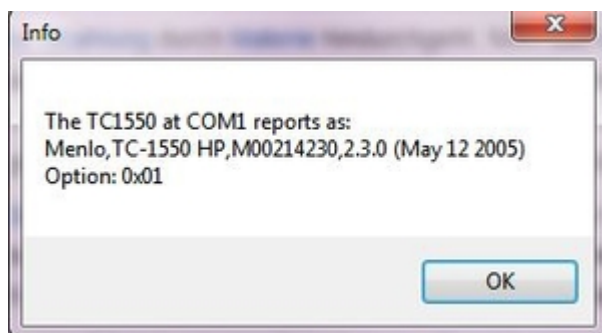


- If you **turn off and on** the frequency comb, then push in the order:
 1. *Reset*
 2. *Auto on*
 3. *Enable on*



The program is done, if all four boxes [*Laser on*, *Modelocked*, *Autoscrumble*, *Enabled*] are green!

- For **closing** the software:
 1. *Halt*
 2. *Freeze*
 3. Close the window

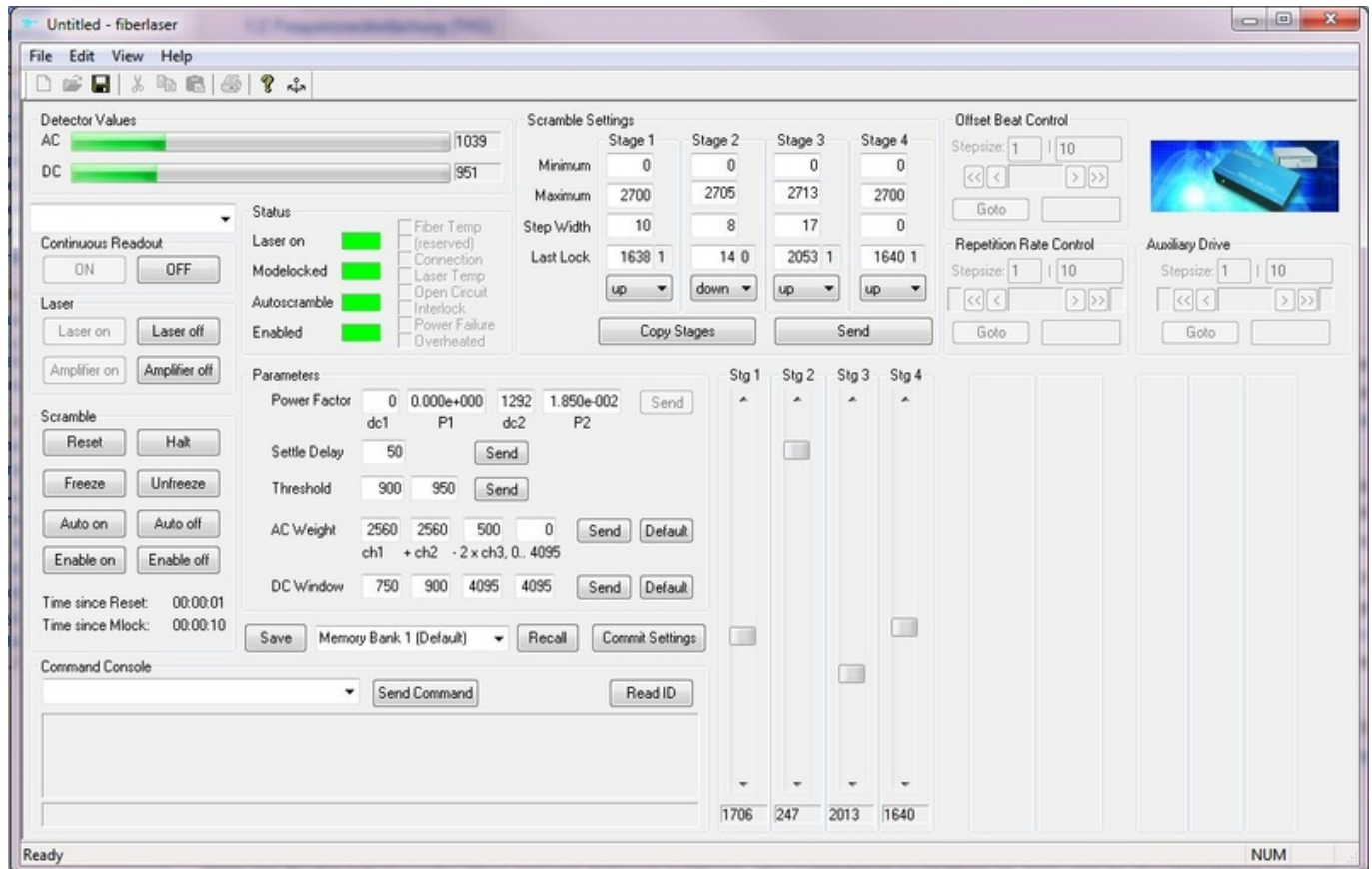

From:

<https://iqwiki.iqo.uni-hannover.de/> - IQwiki

Permanent link:

<https://iqwiki.iqo.uni-hannover.de/doku.php?id=groups:mg:fiberlaser&rev=1481624883>

Last update: **2016/12/13 10:28**

