



If you are present at a power-failure, prepare the Emergency-Power-System (in D320 Mg-Lab next to the electronics/chemistry cabinet) and get gasoline (Super Benzin)

At power failure

- Set up the emergency power system (EPS) on the balcony room - (Parameters to turn it on: Choke: Turn off, Gas: Turn on)
- Disconnect the marked power-cord (next to "Notaus") from the power-plug and put it into the EPS. Now, the turbo-pumps and IGP should be save
- If the turbo-pumps fail: Close the big valve below the chamber-table for the big turbo pump and turn off the big turbo.

After power failure



!Most probably the ACs will fail after a powe failure!

- Turn off every device (Oscilloscopes, commercial systems etc) and all the power supplies in the "Stromschrank" to avoid unecessary heating of the lab
- Check if the AC is working again (call 4440 to tell them or wait for the next day until it's fixed).<hi #ed1c24> Indication for a not working AC</hi>: The red light next to the WC/KRb Office is flashing (but musn't)

From:

<https://iqwiki.iqo.uni-hannover.de/> - IQwiki

Permanent link:

<https://iqwiki.iqo.uni-hannover.de/doku.php?id=groups:mg:powerfailure&rev=1484823956>

Last update: **2017/01/19 11:05**

