


TracerDAQ

DAQ Hardware settings

- Optimum: +/- 2V. If there is more voltage, then are the resolution of the fluctuation bad.
- You can change the time-range in the program *TracerDAQ*

Software settings


- Select *Strip Chart*
- Push *Run*

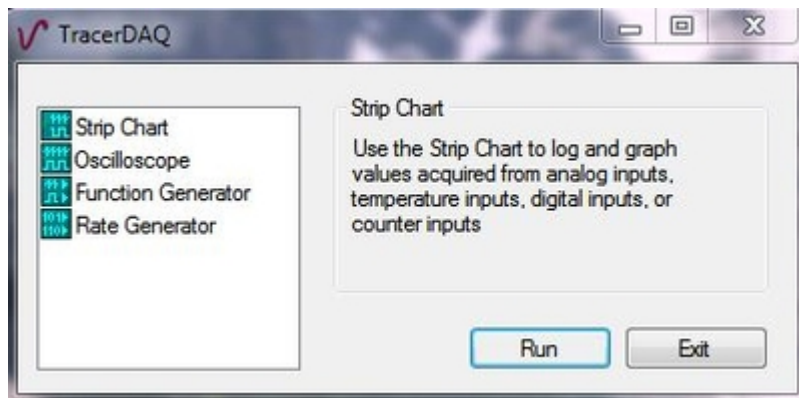


- You can change the range under: Edit → Scan Rate/Trigger Settings
- Push *OK*




The memory will be very fast full for high scan ranges!

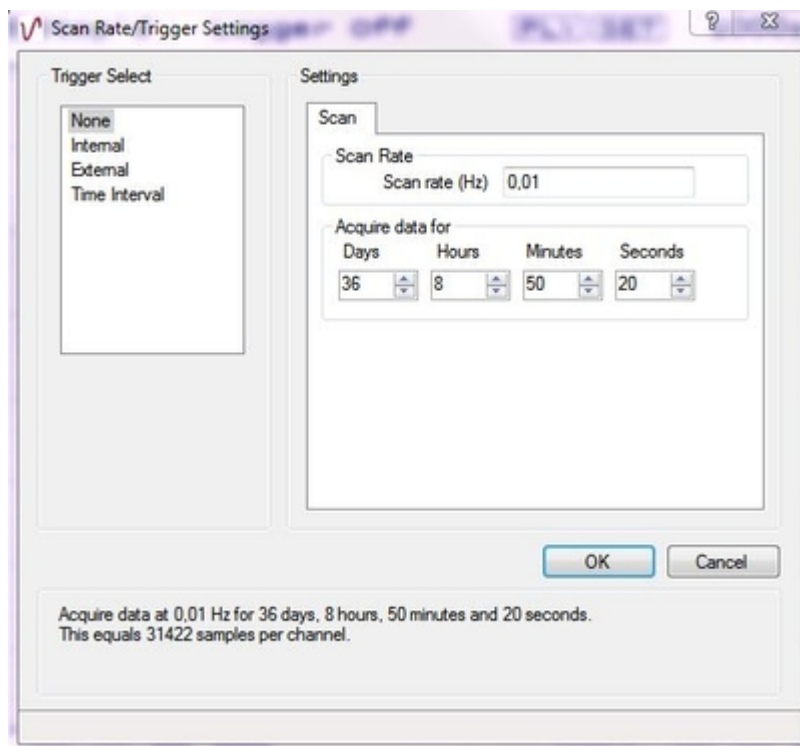

- You have to set the *DAQ Range/Mode* (=Input Voltage) to the right value!
- In this window, you can see the error messages, like *not connected*!

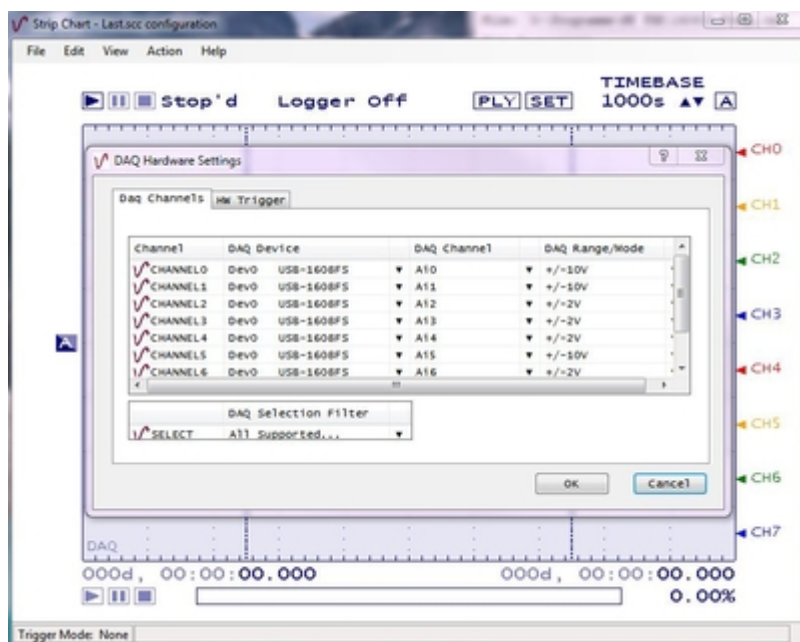

- Push *Play/Start*

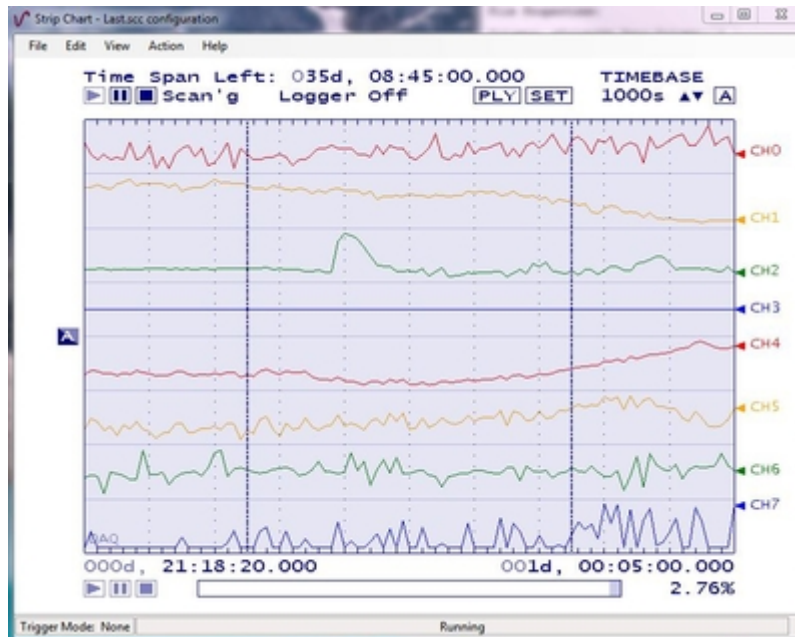

Channels on RedLab-Card:

- 0: R1 Inner T OOL
- 1: R2 Inner T OOL
- 2: R1 PDH Out
- 3: R1 Intens Out
- 4: R2 PDH Out
- 5: R2 Intens Out
- 6: R1 Vacuum
- 7: ECDL

Current Setup: Check Dokumente Folder

From:
<https://iqwiki.iqo.uni-hannover.de/> - IQwiki

Permanent link:
<https://iqwiki.iqo.uni-hannover.de/doku.php?id=groups:mg:private:resonatoren:reslabcomputers:tracerdaq&rev=1578992209>

Last update: 2020/01/14 08:56

